

LED Ring Light (MA341103)

Instruction Manual

Thank you for using our products. This manual is for MA341103 LED Ring Light. Please use the products correctly according to this Instruction Manual.

Specifications:

Power supply: 90-265V 50/60Hz

Output Voltage: 24V

Power Consumption: 6W

Color Temperature: 6400K

Service Life: ≥ 10000 hours

Operating Temperature: 32-104F (0-40°C)

Current: 1A

Max. Opening: 62mm

Quantity of Bulbs: 60 bulbs

Brightness: Adjustable

Operation:

Step 1: Connect to power supply, check if LED Ring Light works normally, and turn it off.

Step 2: Attach the LED light to the microscope.

1. Loosen the three screws on LED
2. Attach the LED light to the microscope
3. Tighten the screws and adjust the LED light

Step 3: Re-start the power supply and adjust the light brightness.



Ring Adapter optional

