

LED Ring Light (MA341305)

Instruction Manual

Thank you for using our products. This Instruction Manual is for MA341305 LED Ring Light. Please use the products correctly according to this Instruction Manual.

Specifications:

Power supply: AC 90-240V 50/60Hz

Max. Opening: 61mm

Quantity of Bulbs: 144 LED Bulbs.

Ring light is divided into 4 areas for quad lighting



Operation

Step 1: Connect to power supply, check if LED Ring Light works normally, and turn it off.



Step 2: Attach the LED light to the microscope.

1. Loosen the three screws on LED
2. Attach the LED light to the microscope
3. Tighten the screws and adjust the LED light



Step 3: According to the different conditions, the ring light is divided into 4 areas and each area is turned on and off separately. Brightness is adjustable.

