

LED Ring Light (MA342103)

Instruction Manual

Thank you for using our products. This Instruction Manual is for MA342103 LED Ring Light. Please use the products correctly according to this Instruction Manual.

Specification:

Power supply: 90-265V 50/60Hz


Quantity of Bulbs: 96 bulbs

Power Output: 4W

I.D.: 50mm, O.D.: 68mm

Brightness: Adjustable

Fits for MZ, MZ-II and MZ-III series microscope bodies.

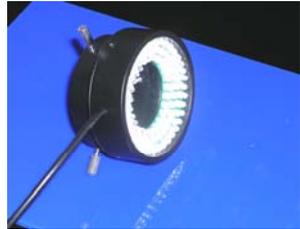

Operation:

Step 1: Connect to power supply, check if LED Ring Light works normally, and turn it off.



Step 2: Attach the LED light to the microscope.

1. Loosen the three screws on LED
2. Attach the LED light to the microscope
3. Tighten the screws and adjust the LED light

Step 3: Re-start the power supply and adjust the light brightness.